

## Product Information

### Caspase-3 Colorimetric Assay Kit

#### I. Kit Contents:

Component	K2008-25 25 assays	K2008-100 100 assays	K2008-200 200 assays	K2008-400 400 assays	Part Number
Cell Lysis Buffer	25 ml	100 ml	100 ml	100 ml	K2008-C-1
2X Reaction Buffer	2 ml	4 x 2 ml	16 ml	32 ml	K2008-C-2
DEVD-pNA (4 mM)	125 µl	500 µl	2 x 0.5 ml	2 x 1 ml	K2008-C-3
DTT (1 M)	100 µl	400 µl	400 µl	400 µl	K2008-C-4
Dilution Buffer	25 ml	100 ml	200 ml	400 ml	K2008-C-5

#### II. Introduction:

Cysteine-dependent aspartate-directed proteases (Caspases) are a family of cysteine proteases that play important roles in apoptosis, necrosis, and inflammation. Sequential activation of caspases plays an important role in cell apoptosis. Caspase-3 is a caspase protein that cleaves and activates Caspases-6 and -7, and is processed and activated by Caspases-8, -9, and 10. Caspase-3 is the predominant caspase involved in the cleavage of amyloid-beta 4A precursor protein, which is associated with Alzheimer's disease. Caspase-3 recognizes tetra-peptide sequences D-x-x-D and hydrolyzes peptide bonds after aspartic acid residues.

Caspase-3 Colorimetric Assay Kit provides a convenient and simple way for detecting the DEVD-dependent caspase activity. When cleavage of DEVD from the labeled substrate DEVD-p-nitroaniline (DEVD-pNA) by Caspase-3 or related caspases, the free pNA light emission can be quantified by using a microtiter plate reader or a spectrophotometer at 405 or 400 nm. Comparison of the absorbance of free pNA from an apoptotic sample with an uninduced control determines the fold increase in Caspase-3 activity.

#### III. Caspase-3 Assay Protocol:

##### A. General Considerations

Aliquot enough 2X Reaction Buffer for the number of assays to be performed. Add DTT to the 2X Reaction Buffer immediately before use (10 mM final concentration: add 10 µl of 1.0 M DTT stock per 1 ml of 2X Reaction Buffer).

Protect DEVD-pNA from light.

##### B. Assay Procedure

1. Induce apoptosis in cells by desired method. Concurrently incubate a control culture without induction.
2. Count cells and pellet  $1-5 \times 10^6$  cells.
3. Resuspend cells in 50 µl of chilled Cell Lysis Buffer and incubate cells on ice for 10 minutes.
4. Centrifuge for 1 min in a microcentrifuge (10,000 x g).
5. Transfer supernatant (cytosolic extract) to a fresh tube and put on ice for immediate assay or aliquot and store at  $-80^{\circ}\text{C}$  for future use.
6. Assay protein concentration.
7. Dilute 50-200 µg protein to 50 µl Cell Lysis Buffer for each assay.
8. Add 50 µl of 2X Reaction Buffer (containing 10 mM DTT) to each sample.
9. Add 5 µl of the 4 mM DEVD-pNA substrate (200 µM final conc.) and incubate at  $37^{\circ}\text{C}$  for 1-2 hour.

10. Read samples at 400- or 405-nm in a microtiter plate reader, or spectrophotometer using a 100- $\mu$ l micro quartz cuvette (Sigma), or dilute sample to 1 ml with Dilution Buffer and using regular cuvette (note: Dilution of the samples proportionally decreases the reading). Fold-increase in CPP32 activity can be determined by comparing these results with the level of the uninduced control.

### Notes:

Background reading from cell lysates and buffers should be subtracted from the readings of both induced and the uninduced samples before calculating fold increase in CPP32 activity.

### General Troubleshooting Guide for Caspase Colorimetric and Fluorometric Kits:

Problems	Cause	Solution
Assay not working	<ul style="list-style-type: none"> <li>• Cells did not lyse completely</li> <li>• Experiment was not performed at optimal time after apoptosis induction</li> <li>• Plate read at incorrect wavelength</li> <li>• Old DTT used</li> </ul>	<ul style="list-style-type: none"> <li>• Resuspend the cell pellet in the lysis buffer and incubate as described in the datasheet</li> <li>• Perform a time-course induction experiment for apoptosis</li> <li>• Check the wavelength listed in the datasheet and the filter settings of the instrument</li> <li>• Always use freshly thawed DTT in the cell lysis buffer</li> </ul>
High Background	<ul style="list-style-type: none"> <li>• Increased amount of cell lysate used</li> <li>• Increased amounts of components added due to incorrect pipetting</li> <li>• Incubation of cell samples for extended periods</li> <li>• Use of expired kit or improperly stored reagents</li> <li>• Contaminated cells</li> </ul>	<ul style="list-style-type: none"> <li>• Refer to datasheet and use the suggested cell number to prepare lysates</li> <li>• Use calibrated pipettes</li> <li>• Refer to datasheet and incubate for exact times</li> <li>• Always check the expiry date and store the individual components appropriately</li> <li>• Check for bacteria/ yeast/ mycoplasma contamination</li> </ul>
Lower signal levels	<ul style="list-style-type: none"> <li>• Cells did not initiate apoptosis</li> <li>• Very few cells used for analysis</li> <li>• Use of samples stored for a long time</li> <li>• Incorrect setting of the equipment used to read samples</li> <li>• Allowing the reagents to sit for extended times on ice</li> </ul>	<ul style="list-style-type: none"> <li>• Determine the time-point for initiation of apoptosis after induction (time-course experiment)</li> <li>• Refer to datasheet for appropriate cell number</li> <li>• Use fresh samples or aliquot and store and use within one month for the assay</li> <li>• Refer to datasheet and use the recommended filter setting</li> <li>• Always thaw and prepare fresh reaction mix before use</li> </ul>
Samples with erratic readings	<ul style="list-style-type: none"> <li>• Uneven number of cells seeded in the wells</li> <li>• Samples prepared in a different buffer</li> <li>• Adherent cells dislodged and lost at the time of experiment</li> <li>• Cell/ tissue samples were not completely homogenized</li> <li>• Samples used after multiple freeze-thaw cycles</li> <li>• Presence of interfering substance in the sample</li> <li>• Use of old or inappropriately stored samples</li> </ul>	<ul style="list-style-type: none"> <li>• Seed only equal number of healthy cells (correct passage number)</li> <li>• Use the cell lysis buffer provided in the kit</li> <li>• Perform experiment gently and in duplicates/triplicates; apoptotic cells may become floaters</li> <li>• Use Dounce homogenizer (increase the number of strokes); observe efficiency of lysis under microscope</li> <li>• Aliquot and freeze samples, if needed to use multiple times</li> <li>• Troubleshoot as needed</li> <li>• Use fresh samples or store at correct temperatures until use</li> </ul>
Unanticipated results	<ul style="list-style-type: none"> <li>• Measured at incorrect wavelength</li> <li>• Cell samples contain interfering substances</li> </ul>	<ul style="list-style-type: none"> <li>• Check the equipment and the filter setting</li> <li>• Troubleshoot if it interferes with the kit (run proper</li> </ul>

		controls)
General issues	<ul style="list-style-type: none"> <li>• Improperly thawed components</li> <li>• Incorrect incubation times or temperatures</li> <li>• Incorrect volumes used</li> <li>• Air bubbles formed in the well/tube</li> <li>• Substituting reagents from older kits/ lots</li> <li>• Use of a different 96-well plate</li> </ul>	<ul style="list-style-type: none"> <li>• Thaw all components completely and mix gently before use</li> <li>• Refer to datasheet &amp; verify the correct incubation times and temperatures</li> <li>• Use calibrated pipettes and aliquot correctly</li> <li>• Pipette gently against the wall of the well/tubes</li> <li>• Use fresh components from the same kit</li> <li>• Fluorescence: Black plates; Absorbance: Clear plates</li> </ul>
<p>Note# The most probable cause is listed under each section. Causes may overlap with other sections.</p>		

**For research use only! Not to be used in humans.**

### **Our promise**

If the product does not perform as described on this datasheet, we will offer a refund or replacement. For more details, please visit <http://www.apexbt.com/> or contact our technical team.

Tel: +1-(832)696-8203

Fax: +1-832-641-3177

Email: [sales@apexbt.com](mailto:sales@apexbt.com)