

Product Information

Fluorescein Active Caspase-8 Staining Kit

I. Kit Contents:

Component	K2050-25	K2050-100	Part Number
	25 assays	100 assays	
FITC-IETD-FMK	25 μ l	100 μ l	K2050-C-1
Wash Buffer	50 ml	2 x 100 ml	K2050-C-2
Z-VAD-FMK	10 μ l	10 μ l	A1902

II. Introduction:

Caspases activation is crucial in apoptosis. Using Fluorescein Caspase-8 Staining Kit is an easy and sensitive way of detecting activated caspase-3 in living cells. This assay uses the Caspase-8 inhibitor, IEVD-FMK, coupled to FITC (FITC-IEVD-FMK) as a marker. In apoptotic cells, the cell permeable and nontoxic FITC-IEVD-FMK irreversibly binds to activated Caspase-8.

III. Caspase-8 Assay Protocol:

A. Staining Procedure:

1. Induce apoptosis in cells (1×10^6 /ml) by desired method. Concurrently incubate a control culture without induction. An additional negative control can be prepared by adding the caspase inhibitor Z-VAD-FMK at 1 μ l/ml to an induced culture to inhibit caspase activation.
2. Aliquot 300 μ l each of the induced and control cultures into eppendorf tubes.
3. Add 1 μ l of FITC-IETD-FMK into each tube and incubate for 0.5-1 hour at 37°C incubator with 5% CO₂.
4. Centrifuge cells at 3000 rpm for 5 minutes and remove supernatant.
5. Resuspend cells in 0.5 ml of Wash Buffer, and centrifuge again.
6. Repeat Step 5.

Proceed to B, C, or D depending on methods of analysis.

B. Quantification by Flow Cytometry:

For flow cytometric analysis, resuspend cells in 300 μ l of Wash buffer. Put samples on ice. Analyze samples by flow cytometry using the FL-1 channel.

C. Detection by Fluorescence Microscopy:

For fluorescence microscopic analysis, resuspend cells in 100 μ l Wash buffer. Add one drop of the cell suspension onto a microslide and cover with a coverslip. Observe cells under a fluorescence microscope using FITC filter. Caspase-8 positive cells appear to have brighter green signals, whereas Caspase-8 negative control cells show much weaker signal.

D. Analysis by Fluorescence Plate Reader: For analysis with fluorescence plate reader, resuspend cells in 100 μ l Wash Buffer and then transfer the cell suspension to each well of the black microtiter plate. Measure the fluorescence intensity at Ex/Em = 485/535 nm. For control, use wells containing unlabeled cells.

General Troubleshooting Guide:

Problems	Cause	Solution
High Background	<ul style="list-style-type: none"> • Cell density is higher than recommended • Cells were not washed well with wash buffer after staining 	<ul style="list-style-type: none"> • Refer to data sheet and use the suggested cell number • Use the wash buffer provided, and as instructed in the

	<ul style="list-style-type: none"> • Cells were Incubated for extended period of time • Use of extremely confluent cells • Contaminated cells 	<p>datasheet</p> <ul style="list-style-type: none"> • Refer to data sheets and incubate for exact times • Perform assay when cells are at 70-95% confluency • Check for bacteria/ yeast/ mycoplasma contamination
Lower signal levels	<ul style="list-style-type: none"> • Cells did not initiate apoptosis • Very few cells used for analysis • Incorrect setting of the equipment used to read samples • Use of expired kit or improperly stored reagents 	<ul style="list-style-type: none"> • Determine the time-point for initiation of apoptosis after induction (time-course experiment) • Refer to data sheet for appropriate cell number • Refer to data sheet and use the recommended filter setting • Always check the expiry date and store the components appropriately
Erratic results	<ul style="list-style-type: none"> • Old (unhealthy) cells used • Adherent cells were dislodged and washed away prior to assaying • Incorrect incubation times or temperatures • Incorrect volumes used 	<ul style="list-style-type: none"> • Seed healthy cells and make sure cells are healthy prior to induction of apoptosis • Collect all cells (both attached and dislodged) after induction for accurate results • Refer to datasheet & verify correct incubation times and temperatures • Use calibrated pipettes and aliquot correctly

Note: The most probable cause is listed under each section. Causes may overlap with other sections

For research use only! Not to be used in humans.

Our promise

If the product does not perform as described on this datasheet, we will offer a refund or replacement. For more details, please visit <http://www.apexbt.com/> or contact our technical team.

Tel: +1-(832)696-8203

Fax: +1-832-641-3177

Email: sales@apexbt.com