

Product Information

Hydrogen Peroxide Assay Kit

I. Kit Contents:

Components	K2110-200 200 assays	Cap Color	Part Number
H ₂ O ₂ Assay Buffer	25 ml	WM	K2110-C-1
OxiRed™ Probe (in anhydrous DMSO)	0.2 ml	Red	K2110-C-2
HRP	1 vial	Green	K2110-C-3
H ₂ O ₂ Standard (0.88 M)	0.1 mg	Yellow	K2110-C-4

II. Introduction:

Hydrogen Peroxide plays an important role in various oxidative stress-associated conditions. Hydroperoxide-mediated NF-κB is related to asthma, neurodegenerative disease, atherosclerosis and immune system etc. Hydrogen Peroxide Assay Kit provides a sensitive, easy and direct way for measuring H₂O₂ in biological samples. In the presence of Horse Radish Peroxidase (HRP), the OxiRed Probe reacts with H₂O₂ to produce product with color (λ_{max} = 570 nm) and red-fluorescent (Ex/Em = 535/587 nm). The detection limit can be as low as 2 pmol per assay (or 40 nM concentration) of H₂O₂ in the sensitive fluorometric assay.

III. Storage and Handling:

Warm the assay buffer to room temperature before use. Briefly centrifuge vials prior to opening. Read the entire protocol before performing the assay.

IV. Reagent Reconstitution and General Consideration:

OxiRed™ Probe: Ready to use as supplied. Briefly warm at 37°C to melt frozen DMSO. The OxiRed™ Probe solution is stable for 1 week at 4°C and 1 month at -20°C.

HRP: Dissolve in 220 µl Assay Buffer, pipetting up and down. The HRP solution is stable for 1 week at 4°C and 1 month at -20°C.

V. Hydrogen Peroxide Assay Protocol:

- Sample Preparations:** Collect cell culture supernatant, serum, plasma, urine and other biological fluids (contains 0.8 - 6 µM H₂O₂). Centrifuge for 15 minutes at 1000 x g within 30 minutes of collection. Remove particulate pellet. Samples, especially those such as culture medium, tissue lysate or plasma should be filtered through a 10 kDa MW spin filter to remove all proteins then kept at -80°C for storage. It is recommended with all sample types to assay immediately or aliquot and store the samples at -80°C. Avoid repeated freeze-thaw cycles. Add 2 - 50 µl samples into each well, bring the volume to 50 µl with assay buffer.
- H₂O₂ Standard Curve:** For the Colorimetric Assay: Dilute 10 µl 0.88 M H₂O₂ standard into 870 µl dH₂O to generate a 10 mM H₂O₂ standard, then dilute 10 µl of the 10 mM H₂O₂ standard into 990 µl dH₂O to generate a 0.1 mM H₂O₂ standard. Add 0, 10, 20, 30, 40, 50 µl of the 0.1 mM H₂O₂ standard into 96-well plate in duplicate to generate 0, 1, 2, 3, 4, 5 nmol/well H₂O₂ standard. For the Fluorometric Assay: Dilute 100 µl of the 0.1 mM

H₂O₂ standard into 900 µl dH₂O to generate a 10 µM H₂O₂ Standard. Add 0, 10, 20, 30, 40, 50 µl of the 10 µM H₂O₂ standard into a 96-well plate in duplicate to generate 0, 0.1, 0.2, 0.3, 0.4, 0.5 nmol/well H₂O₂ standard.

3. Reaction Mix: Mix enough reagents for the number of assays to be performed. For each well, prepare a total 50 µl Reaction Mix:

	Colorimetric Assay	Fluorometric Assay
Assay Buffer	46 µl	48 µl
OxiRed™ Probe solution	2 µl	1 µl
HRP solution	2 µl	1 µl

Add 50 µl of the Reaction Mix to each test samples and H₂O₂ standards. Mix well. Incubate at room temperature for 10 min.

For a more sensitive assay, you can dilute the standard 10 fold further, decrease OxiRed™ amount to 0.2 µl and HRP amount to 0.4 µl per well, it will decrease the fluorescence background and detects as low as 2 pmol/well (or 40 µM) H₂O₂.


4. Measure OD (570 nm) or fluorescence (Ex/Em = 535/587 nm) in a micro-plate reader.

5. Calculation: Correct background by subtracting the value derived from the 0 nmol H₂O₂ control from all sample and standard readings (Note: The background reading can be significant and must be subtracted from sample readings). Plot the H₂O₂ standard curve. Apply your sample readings to the standard curve. H₂O₂ concentrations of the test samples can then be calculated,

$$C = Sa/Sv \text{ (pmol/}\mu\text{l or }\mu\text{M)},$$

Where: Sa is the sample amount from your standard curve (in pmol),

Sv is sample volume (µl).

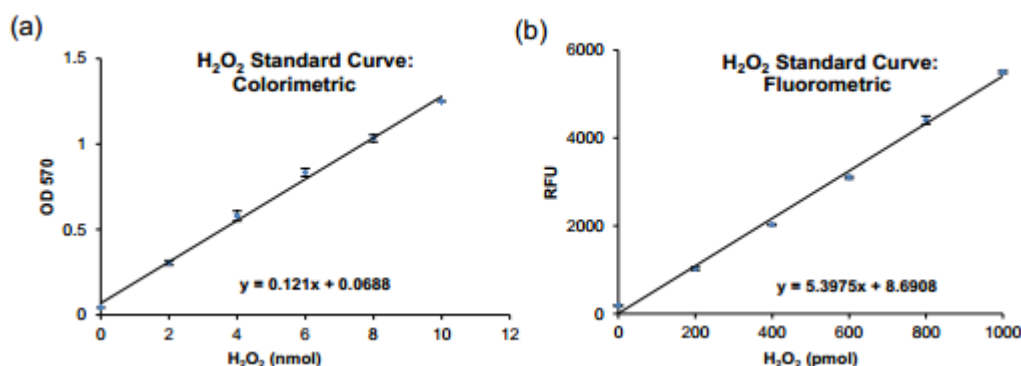

General Troubleshooting Guide:

Problems	Cause	Solution
Assay not working	<ul style="list-style-type: none"> • Use of ice-cold assay buffer • Omission of a step in the protocol • Plate read at incorrect wavelength • Use of a different 96-well plate 	<ul style="list-style-type: none"> • Assay buffer must be at room temperature • Refer and follow the data sheet precisely • Check the wavelength in the data sheet and the filter settings of the instrument • Fluorescence: Black plates ; Luminescence: White plates; Colorimeters: Clear plates
Samples with erratic readings	<ul style="list-style-type: none"> • Use of an incompatible sample type • Samples prepared in a different buffer • Samples were not deproteinized (if indicated in datasheet) • Cell/ tissue samples were not completely homogenized • Samples used after multiple free-thaw cycles • Presence of interfering substance in the sample 	<ul style="list-style-type: none"> • Refer data sheet for details about incompatible samples • Use the reaction buffer provided in the kit or refer data sheet for instructions • Use the 10 kDa spin cut-off filter or PCA precipitation as indicated • Use Dounce homogenizer (increase the number of strokes);

	<ul style="list-style-type: none"> • Use of old or inappropriately stored samples 	observe for lysis under microscope <ul style="list-style-type: none"> • Aliquot and freeze samples if needed to use multiple times • Troubleshoot if needed, deproteinize samples • Use fresh samples or store at correct temperatures till use
Lower/ Higher readings in Samples and Standards	<ul style="list-style-type: none"> • Improperly thawed components • Use of expired kit or improperly stored reagents • Allowing the reagents to sit for extended times on ice • Incorrect incubation times or temperatures • Incorrect volumes used 	<ul style="list-style-type: none"> • Thaw all components completely and mix gently before use • Always check the expiry date and store the components appropriately • Always thaw and prepare fresh reaction mix before use • Refer data sheet & verify correct incubation times and temperatures • Use calibrated pipettes and aliquot correctly
Readings do not follow a linear pattern for Standard curve	<ul style="list-style-type: none"> • Use of partially thawed components • Pipetting errors in the standard • Pipetting errors in the reaction mix • Air bubbles formed in well • Standard stock is at an incorrect concentration • Calculation errors • Substituting reagents from older kits/ lots 	<ul style="list-style-type: none"> • Thaw and re-suspend all components before preparing the reaction mix • Avoid pipetting small volumes • Pipette gently against the wall of the tubes • Prepare a master reaction mix whenever possible • Always refer the dilutions in the data sheet • Recheck calculations after referring the data sheet • Use fresh components from the same kit
Unanticipated results	<ul style="list-style-type: none"> • Measured at incorrect wavelength • Samples contain interfering substances • Use of incompatible sample type • Sample readings above/below the linear range 	<ul style="list-style-type: none"> • Check the equipment and the filter setting • Troubleshoot if it interferes with the kit • Refer data sheet to check if sample is compatible with the kit or optimization is needed • Concentrate/ Dilute sample so as to be in the linear range
Note: The most probable list of causes is under each problem section. Causes/ Solutions may overlap with other problems.		

For research use only! Not to be used in humans.

Our promise

If the product does not perform as described on this datasheet, we will offer a refund or replacement. For more details, please visit <http://www.apexbt.com/> or contact our technical team.

Tel: +1-(832)696-8203
 Fax: +1-832-641-3177
 Email: sales@apexbt.com