

Protocol Cat. No. K3012

## **BstBl**

### Introduction

BstBl is a genetically engineered fast restriction endonuclease suitable for fast digestion of plasmid DNA, PCR products or genomic DNA, etc. (enzyme digestion is completed within 5-15 minutes). Its internal matching Reaction Buffer (Color) contains red and yellow tracer dyes for direct use in gel electrophoresis. The migration rate of red dye and 2500 bp double-stranded DNA fragment in 1% agarose gel was similar, and the migration rate of yellow dye and 10 bp double-stranded DNA fragment in 1% agarose gel was similar.

Its recognition sequence and cutting site are as follows:

5'.....TT<sup>1</sup>CGAA.....3'

3'......5'

# Components and Storage

	Unoun
Espore I	e Un
Peresterion.	
a service and the service and	

Size	K3012 - 2000 U	Storage	
Components			
BstBI (20 U/µL)	100 µL	-20°C	
10×Reaction Buffer	1 mL	-20°C	
10×Reaction Buffer (Color)	1 mL	-20°C	
Shipping: Dry Ice	Shelf life: 2 years	E good mound	
Dratter of the second		Perfection	

## Protocol

#### I DNA Rapid enzyme digestion:

1. Refer to the table below to configure the reaction system on the ice:

Components	Plasmid DNA	PCR products	Genomic DNA
DNA substrate	Х µL (1 µg)	X μL (~0.2 μg)	Х µL (5 µg)
10×Reaction Buffer or		2.11	Expore the On
10×Reaction Buffer (Color)	μ Ζμι		erection 5 µL
BstBl	1 µL	1 μL	5 µL
Nuclease-free Water	Το 20 μL	Το 30 μL	To 50 μL
Total Reaction Volume	20 µL	30 µL	50 µL

**Note:** This system is suitable for enzymatic digestion of purified PCR products; Since DNase also has exonuclidene activity, it will affect the enzyme digestion products, so the following step needs to be cloned and other operations, it is recommended to purify the PCR products before enzyme digestion.

- 2. Gently mix the reaction system (gently suck or flick the tube wall, do not swirl), and then instantaneously centrifuge to collect the tube wall residue.
- 3. Incubate at 37 °C for 15 min (plasmid); 15-30 min (PCR product); 30-60 min (genomic DNA).
- Heating at 80°C for 20 min inactivates BstBI and terminates the reaction (optional). 4.
- 5. If 10×Reaction Buffer (Color) is selected for enzyme digestion, the obtained product can be directly subjected to sample electrophoresis.

#### Il Double or multiple enzyme digestion:

- 1. The dosage of each endonuclide is 1 µL, and the reaction system is appropriately expanded as needed.
- 2. The total volume of all endonuclease enzymes should not exceed 1/10 of the total reaction system.

3. If the optimal reaction temperature of several selected endonucleoenzymes is different, the enzyme with the lowest optimal temperature should be started first, and then the optimal temperature should be added

PEX

For higher enzymes, incubate at higher temperatures.

#### **Notes**

- 1, Add the BstBI enzyme last when configuring the enzyme digestion reaction system, and keep the BstBI on ice when removing it from the refrigerator.
- If the total reaction system is greater than 20 µL, the incubation time should be appropriately increased, and 2. the water bath, metal bath or sand bath should be used as far as possible.
- Mix the components by blowing up and down the pipette or "flicking" the reaction tube, do not use violent 3. vortex mixing.
- 4. For DNA methylated by CpG or EcoKI, clipping may be blocked; In addition, asterisk activity may occur over a incubation time of 3 h.
- 5. For double or multi-enzyme digestion, an appropriate buffer compatible with two or more endonuclide enzymes should be selected, and the reaction system should be set up according to the table above. If there is no suitable buffer to choose from, one enzyme can be digested and purified first, and then another enzyme APEABLE BIL digestion reaction can be performed after purification.
- 6. For research use only. Not to be used in clinical diagnostic or clinical trials.

