

# HyperTrap Phenyl HP Column

## Product description

HyperChrom Phenyl HP Agarose is a class of hydrophobic interaction chromatography media that separates molecules using differences in their hydrophobic properties and how much they are under specific conditions. The HyperChrom Phenyl HP Agarose series has excellent scale-up performance: 1) Highly crosslinked agarose base frame with excellent rigidity, so high process flow rates and improved process efficiency can be achieved at low backpressure. 2) Hydrophilic base frame to minimize the influence on the hydrophobic role of the ligand. 3) Through chemical modification, it has excellent chemical compatibility and is resistant to CIP cleaning such as sodium hydroxide. 4) Fine particle size design to improve resolution.

HyperTrap Phenyl HP Column is a preloaded column with HyperChrom Phenyl HP Agoros as the chromatography medium, which is mainly used for the isolation and purification of a variety of biomolecules, such as: recombinant proteins, antibodies, viral vaccines, etc.

The main body and inner plug of the pre-installed column tube are made of PP, and the inside and outside are polished, and the sieve plate is made of HDPE material, which has excellent physical properties and chemical resistance, corrosion resistance, anti-aging, long life and good safety. 1) Ready-to-use, convenient and fast. 2) Can be used with syringes, peristaltic pumps or chromatography systems. 3) It can be used in series to increase the sample processing capacity.

# APENER BIO

## Product parameters

Pre-assembled column specification/name	1 mL	5 mL		
Dimensions (IDxH)	7.7 x 25 mm	16 x 25 mm		
Withstand pressure	0.3 MPa	0.3 MPa		
Workflow Speed*	1 mL/min	1-3 mL/min		
Operating temperature	4 - 30°C	4 - 30°C		
Storage condition	Store the components at 4°C for 5 years.			
*This flow rate is the recommended common flow rate, the pressure resistance, and the use of different				
types of chromatography packing in the pre-assembled column will be different, please refer to the				
product manual of the corresponding chromatographic filler.				

1. HyperTrap Phenyl HP Column preloaded column parameters

2. HyperTrap Phenyl HP Column chromatography media parameters

Name	Description	
Chromatograp	Hydrophobic interactions	
hy media type		
Ligation	R-O-CH2-CH(OH)-CH2 -O-C6H5	
Scaffolding	Highly cross-linked agarose	
Average	34 um	
particle size		
Ligand density	25 μmol phenyl group/mL chromatography medium	
Dynamic load	>45 mg α-chymotrypsinogen/mL chromatography medium*	
Flow rates are	90-150 cm/h	
recommended		
Maximum	200 cm/h	
flow rate		
Withstand	0.3 MPa	
pressure		
Use	1 20%	
temperature	4-30 C	
pH stability**	2-14 grander	
Solvent resistant	Common aqueous solution, 3 M ammonium sulfate, 30% isopropanol***, 75% ethanol***, 10% ethylene glycol***, 1 M NaOH, 1 M acetic acid, 6 M guanidine hydrochloride, 8 M urea	
Intolerant to solvents	Oxidant	

\*Dynamic load measurement conditions: column height: 10 cm, test flow rate 150 cm/h, test buffer: 0.1 M potassium phosphate, 2 M ammonium sulfate solution, pH 7.0, test sample: 2mg/m L of  $\alpha$ -chymotrypsinogen sample, when the penetration of  $\alpha$ -chymotrypsinogen reaches 10%, the amount of  $\alpha$ -chymotrypsinogen per unit volume (mL) (mg).

\*\*After 7 days in the environment of chromatography at 40°C and pH 2-14, the physical and chemical properties and functions of the chromatography medium did not change significantly.

\*\*\* 30%, 75% and 10% are v/v, volume ratio.

# Experimental manipulation

#### 1. Preparation of buffers

Buffer selection: Buffer salts that do not interact with the chromatography medium should be selected.

- If the binding elution mode is used: a high pH buffer (usually 1 pH unit higher than the isoelectric point of the target molecule) buffer should be used for the equilibrium buffer to facilitate the binding of the target molecule, and the stability of the sample in the buffer should be considered; Elution buffers are usually buffers in which a high concentration of salt (e.g., 1 M NaCl) is added to the equilibration buffer.
- If the flow through mode is used: the equilibration buffer should be treated under conditions conducive to the binding of impurities, and rinsed directly with a high concentration of salt after the target molecule has completely flowed through.

#### 2. Sample preparation

Prepare samples for purification.

#### 3. Chromatographic conditions

- Flow rate selection: Linear flow rate of 90-150 cm/h is generally selected according to the height of the column bed.
- Sample preparation: To prevent the sample from clogging the column, the sample needs to be filtered with a 0.2/0.45  $\mu$ m (after inclusion body disruption) microporous membrane before loading, and it is recommended that the pH and conductivity of the sample be adjusted to be consistent with the equilibrium buffer (the pH and conductivity of the sample can be adjusted by dilution, ultrafiltration, and desalting)  $_{\circ}$

#### 4. Preloaded column setup

- (1) Open the package and remove the preloaded column.
- (2) Connect the pre-packed column: unscrew the upper and lower plugs of the pre-packed column and connect it to the chromatography system (peristaltic pump or syringe, etc.), rinse the pre-packed column with pure water 2-3 times the column volume to drain the preservation solution (generally 20% ethanol). In order to prevent air bubbles from entering, pure water can be dripped and filled at the inlet end of the pre-loaded column before connecting the system connector.
- (3) Cleaning and disinfection: For the first use, it is recommended to clean and disinfect the column and rinse 2 times the column volume with pure water or buffer. For recommended cleaning and disinfection buffers, refer to subsequent cleaning and regeneration procedures.
- (4) Sample preparation: To prevent the sample from clogging the column, it is recommended that the sample be filtered with a 0.45 μm microporous membrane before loading.
- 5. Column efficiency determination (optional).

Select one of the two test methods shown in the table below for column effectiveness testing. Use the mobile phase equilibrium chromatography column to the baseline to be stable, load the sample into the chromatography column, continue to use the mobile phase for rinsing, and after the chromatographic peak is completed to return to the baseline, end the run, integrate the chromatographic peak, and evaluate the loading effect.

	Acetone method	NaCl method
Sample	1% (v/v) acetone in water	2 M NaCl in water
Sample volume	1% column volume	1% column volume
Mobile phase	water	0.2 M NaCl in water

Table 2 Statistical table of two column efficiency measurement methods

Velocity of flow	30 cm/h	30 cm/h
Detector	UV 280 nm	electrical conductivity

The main evaluation criteria for the effect of column loading are N/m (number of plates per meter) and As (symmetry factor), which are calculated as follows:



CAPENER CONT

Column efficiency qualification standards: N/m > 3000, 0.8 < As < 1.5

\*Parameter Notes:

L = column height, VR = reserved volume,  $W_h =$  half-peak width, a = left half-peak width at 10% peak height, b = right half-peak width at 10% peak height

#### 6. Chromatographic steps

- Equilibrium: Adequately equilibrate the column to pH and conductivity stable using equilibration buffer and is essentially the same as the equilibration buffer, which typically requires 3-5 times the column volume.
- (2) Sample loading<sup>\*</sup>: Determine the sample loading volume and amount on HyperTrap Phenyl HP Column based on the binding load measured in the pilot experiment.
- (3) Washing<sup>\*</sup>: Rinse the column with equilibration buffer or other suitable buffer until UV stable and return to baseline.
- (4) Elution \*: Elution is achieved by reducing the concentration of salt ions, which can gradually reduce the concentration of salt ions in the eluate through linear gradients or step gradients, and elute molecules with different binding strengths. Fractional collection is performed on the eluted sample. pH gradient elution or mixed elution is also available.
- (5) Regeneration: Rinse the chromatography column with buffer containing low salt.
- (6) Re-equilibration: Re-equilibrate the chromatography column with equilibration buffer.

Note: \* If the flow-through mode is used, the "sample loading" step should be set up for collection; The "washing" step ensures that after all the target molecules have flowed through, the collection can be stopped; In the "elute" step, the impurities are directly eluted with a high-salt buffer.

#### 7. Cleaning and recycling

As the number of uses of the chromatography medium increases, contaminants (e.g., lipids, endotoxins, proteins, etc.) accumulate on the chromatography column. Regular in-place cleaning is essential to keep the column in stable working condition. Determine the frequency of in-place cleaning according to the degree of contamination of the chromatography medium (if the contamination is serious, it is

recommended that in-place cleaning should be carried out after each use to ensure repeatable results and extend the working life of the chromatography medium).

For different types of impurities and contaminants, cleaning can be carried out under the following conditions:

- Removal of strongly binding proteins: Wash with 2 M NaCl solution in 5x column volumes, or use a high salt buffer at pH 2 or higher, such as 1 M NaAc solution.
- Removal of strong hydrophobic proteins and precipitated proteins: first wash with 1 M NaOH solution in 5 column volumes, and then wash the lye with 5-10 column volumes of 20% ethanol.
- Removal of lipoproteins and lipids: Wash first with 5x column volume of 70% ethanol or 30% isopropanol, then rinse with 5-10x column volume of 20% ethanol.

Note: 70% ethanol or 30% isopropanol should be degassed before use; The flow rate during the cleaning process can be selected from 30-60 cm/h; When the blockage is severe, reverse cleaning can be used.

#### 8. Sterilization

In order to reduce the microbial load, it is recommended to use 0.5~1 M NaOH solution to treat the chromatography medium with a processing time of 15~30 min.

#### 9. Stockpile

Unopened chromatography media, please keep in the original container; The completed chromatography column is first soaked with 20% ethanol solution and then the upper and lower column heads are closed. The storage environment is 4~30 °C.

#### **10.** Destruction and recycling

- Since HyperTrap Phenyl HP Column chromatography media is difficult to degrade in nature, incineration of discarded chromatography media is recommended in order to protect the environment.
- For chromatography media exposed to bioactive samples such as viruses and blood, please follow local biosafety requirements before destroying or disposing of them.

#### Notes

- It is recommended that the buffer and protein solution used for purification be filtered through a 0.22 μm or 0.45 μm membrane and then used on the column.
- 2. All chromatography media are kept in 20% ethanol solution with a glue suspension ratio of approximately 75%.
- **3.** This product is for scientific purposes only.

